

Filtration Membranes Must Be Transferred Within The Red Target Area Lines

INCORRECT ❌



CORRECT ✅



Tip: The red lines on the cassette are the "Target Area."

Impact of Error: Incorrect transfer of the membrane to the cassette will result in a Vision Analysis Error (VAE)

Samples Must Be Ordered Prior To Loading On To The System

INCORRECT ❌

Tests	Lot/Batch	Sample ID	Count	Status	Location	Next Image

CORRECT ✅

Tests	Lot/Batch	Sample ID	Count	Status	Location	Next Image
	30oct2019	Dilution2	0	Ordered	UNKNOWN	13:01
	30oct2019	Dilution10	0	Ordered	UNKNOWN	13:01
	30oct2019	Alert	0	Ordered	UNKNOWN	13:01

Tip: Ordered samples will appear on the Home Screen and a Sample ID Label will not be printed.

Impact of Error: Failing to order a sample will prevent the system from knowing what to do with it.

Ensure The Complete 2D Barcode is Printed On The Sample ID Label

INCORRECT ❌



CORRECT ✅



Tip: Use the printer's settings to re-align label stock as necessary. Contact RMB Support for assistance.

Impact of Error: Incomplete barcode printing will prevent the system from identifying a sample and its test parameters.

Apply Sample ID Labels In Designated Area On The Back Of The Cassette

INCORRECT ❌



CORRECT ✅



Tip: Sample ID Labels will fit in molded area on back of the cassette.

Impact of Error: Incorrect label placement will prevent the system from properly reading the barcode to identify the sample's test parameters.

Ensure Vision Lids Are Engaged Correctly

INCORRECT ❌



CORRECT ✅



Tip: The red box on the left shows the locking component of the Vision Lid called a "J-hook" and should not be visible.

Impact of Error: A "Height Check Error" may occur on the system. A Vision Lid may fall off somewhere in the system, requiring service intervention.

Cassettes Must Be Loaded With The Vision Lids Facing Upward

INCORRECT ❌



CORRECT ✅



Impact of Error: The system robotics will be unable to transfer the cassettes within the system.

Make Sure Cassettes Are Loaded Flat In The Carousel



Tip: Ensure that cassettes are stacked flat on top of each other in the carousel. Each carousel has 6 columns that can fit a maximum of 10 cassettes.

Impact of Error: The system robotics will be unable to transfer the cassettes within the system.

Lock The Carousel When Placing Into The System



Tip: Lock the carousel in place by turning the handle clockwise (to the right). You will hear a "click" when the carousel has been locked into place correctly.

Impact of Error: The system robotics will be unable to transfer the cassettes within the system.

All Doors Must Be Opened and Closed When Commanded



System Status

Left Carousel is unlocked

12/20/2019 12:35:34

Growth Direct doors must be opened and closed when commanded via the Home Screen even if you determine you no longer need to open the door after issuing the command.