

Introduction:


This procedure will help you start the Growth Direct® System (GD2). The full referenced document is TSC0010.

Starting the GD System:



Note: These instructions support systems compliant with IEC61010 Version 3. Previously released Growth Direct® Systems have power shutdown steps documented in their user guides.

RMB personnel will install and start the system for you. It is designed to be on all the time. If you lose power, the system will attempt to automatically start when power is restored. However, if you need to manually start the system, follow these steps.

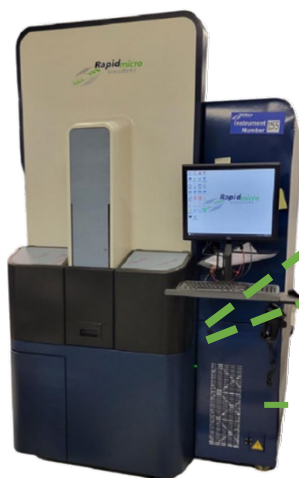

1. Unlock the panel on the bottom right by inserting the panel key (inset) into the lock and turning the key clockwise. The panel swings out exposing the circuit panel.




Lock



Panel Key

Panel

2. If necessary, flip the Main Breaker Switch and UPS Power Switch to their ON positions. These switches are located on the lower right back side of the GD.

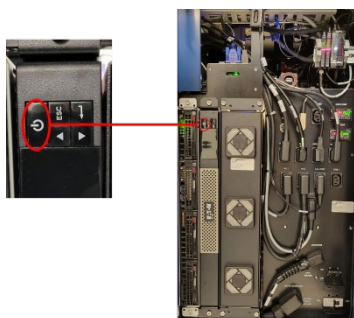

3. Press the UPS Power Button, in the upper left corner of the UPS. All systems, including the servers, monitor, and software turn on automatically.

RMB strives to provide the highest quality products and services to our customers. If you have any questions regarding this informational document, please contact Customer Service at support@rapidmicrobio.com. If interested in placing an order with Rapid Micro Biosystems®, please reach out to orders@rapidmicrobio.com.